Why Use Controlled Nucleation?

Saves money, time and increases product quality. Controlling nucleation during the freezing step in a product to be yophilized has been shown to shorten drying times, decrease reconstitution time and improve the quality traits of the product..

The Controlled Nucleation Advantage

Controlled nucleation of the product eliminates the variation in the degree of super-cooling and consequent freezing rate thereby creating ice crystal growth of optimal pore size to support the fastest sublimation rate possible. Controlled nucleation is designed to create ice matrix uniformity by controlling the freezing of the product in the same way throughout the batch and from cycle to cycle.

Freezing - What Happens if You Don't Have Controlled Nucleation?

The importance of the freezing step is well documented and continues to be studied in the attempt to enhance the characteristics of the frozen matrix prior to lyophilization. Historically the product is "ramped" down in temperature dynamically to a temperature below its critical melting point. Because of a number of factors this simple ramping down in temperature of the product usually creates a great variation in the ice matrix. Most notably the degree of super-cooling is often inconsistent and the size of the resulting ice crystals, in turn, is inconsistent. The ice matrix created during the freezing step essentially creates the resulting pore structure and size, and hence ability of water vapor to leave the product during primary drying. An inconsistent structure will result in inconsistent sublimation rates and cycles must be built conservatively with the worse ice matrix in mind.

How Do I Get Controlled Nucleation on My Existing Laboratory Freeze Dryer?

Just change the door and attach the NS20.

The Millrock FREEZEBOOSTER® is a stand-alone module that can be attached to almost any laboratory freeze dryer. It does not require that the dryer have ASME rated chambers for high pressure and works within the parameters that the normal laboratory lyophilizer is designed for.

FreezeBooster® Controlled Nucleation

Take ultimate control over your freezing process..



FreezeBooster NS20 (portable) attached to a Revo® Lab

Easily retrofit FreezeBooster® on your existing lab freeze dryer.

Will work on lab systems up to 20 sq ft. Installation is as simple as a door swap out FreezeBooster® can be used on mulitiple lyophilizers can be serviced by one FreezeBooster is portable equipment that can be moved from one machine to another.

Avoid hassles, complex and costly factory or field service installation.

No control system upgrades required No additional isolation valves required No external condenser required No ASME rated vessels required

FreezeBooster® Controlled Nucleation

MILLROCK TECHNOLOGY

millrocktech.cor

Millrock's FreezeBooster Controlled Nucleation Station injects ice crystals under the optimum condition into the product chamber to initiate nucleation throughout the entire batch.

FreezeBooster® Controlled Nucleation

FreezeBooster NS20

FEATURES & BENEFITS

On demand nucleation.

Nucleates all vials across the batch at the same time and temperature, resulting in a uniform starting point for crystal growth.

Easily install on any freeze dryer.

One nucleation station can be used on all your lyophilizers with a simple door change. The nucleation station can be used on any freeze dryer brand, and is designed to be moved easily by one operator from lyophilizer to lyophilizer.

Utilize on a varitey of product containers.

The FreezeBooster process has been demonstrated to work on all types of containers, including tubes, microtiter plates, and much more.

Enhance quality & repeatability.

Creates freezing uniformity across the shelf and allows control from batch to batch.

Reduce operating costs.

Lower product cake resistance for shorter primary drying times.

Sterilizable.

Can be sterilized via H2O2, for GMP applications.

PC/PLC controlled system.

System is programmed via a PLC and has the ability for remote access for troubleshooting.

Cost effective.

No need to buy a new freeze dryer No need for major retrofits or system upgrades. Stands alone—one nucleation station will nucleate multiple lyophilizers.

Models



NS20

For use on laboratory style freeze dryers up to 20 sq ft. Interfaces with the product chamber door. Mounted on an adjustable height table, the NS20 version is easily moved between freeze dryers. Option for H2O2 sterilization capability.

NS100

For use with production style lyophilizers rom 30 sq

ft to 100 sq ft. Interfaces with the product chamber door. Console style cabinet that is easily moved between freeze dryers. Option for H2O2 sterilization capability.



NSS100

For use with ASME rated production style freeze dryers from 30 sq ft to 100 sq ft. Interfaces with the product chamber via an 8" port. Skid style format for connection in the service alley. Steam sterilizable.



We also make custom designed units for your larger freeze dryers.

Millrock Technology, Inc 39 Kieffer Lane, Kingston, NY 12401 | 1.845.339.5700 | sales@millrocktech.com Specifications subject to change without noticifation. All specifications based on 20C ambient and 60 Hz Trademarks are registered by Millrock Technology, Inc.



Patent # 8839528, 8875413